Buchi applications Meat & Meat Products





Buchi offers a broad range of NIR applications dedicated to specific segments. The development of a calibration is a matter of experience that includes not only the spectroscopic model itself but the reference analytics in the laboratory. With more than 70 years of practical knowledge Buchi is a competent partner to support your daily business.

NIR Measurements

Measurements can be performed using the NIRFlex N-500 equipped with a Solids Measurement Cell or a Buchi NIRMaster. Standard Petri Dishes are sufficient but a High Performance- or an Unbreakable- Sample Cup will be of advantage in several cases where a maximum of precision and accuracy is requested.

NIR Calibrations

A reliable NIR calibration has to include all kind of variation that is necessarily incorporated into your samples. Our databases consist of thousands of spectra covering geographical and seasonal variation.

Reference Methods

Buchi has many of the laboratory products necessary for an official and accredited analysis of the NIR reference parameters: e.g. heating oven for moisture determination in solids, Soxhlet and Extraction for fat and Kjeldahl for protein content.

Service and Support

NIR spectroscopy is a lively and a living business. Calibration models have to be customized to give the best possible results. Our application specialists will give you the support that you request no matter if you want to use a "Plug and Play" solution with one of our Precalibrations or you want to develop your own chemometric models with our NIRCal software.

Product	Meat	Sausage	Turkey	Chicken Meat	Ham	Pepperoni dryed
	N555-501	N555-502	N555-516	N555-517		uryeu
Property						
Moisture	25.0 - 78.3	22.2 - 80.7	67.0 - 75.0	58.1 - 77.1	52.5 – 64.9	29.7 - 43.2
Protein	5.6 – 26.3	3.9 – 37.1	17.8 - 26.3	14.0 - 24.3	24.1 – 33.1	18.7 - 23.2
Fat	0.8 – 64.0	0.4 – 54.0	1.6 - 15.3	1.3 - 30.5		27.8 - 43.5
Ash		0.6 – 7.5				
Collagen	In development	0.3 – 8.8				
Protein-Collagen		6.3 – 24.9				
Salt					4.2 – 8.8	3.6 - 4.4
Proteolysis Index					19.6 – 36.9	
N (TCA)					5.3 – 11.7	

Parameters:

If not written differently all units are given as [%].

The parameters Proteolysis Index is defined by the ratio: soluble protein / total protein.

N (TCA) refers to the amount of soluble nitrogen determined with Trichloroacetic acid.

Color Codes:

Blue: Buchi Precalibrations

Yellow: Applications in a growing state. Full access to all spectral data with the possibility

for further development

Sample Preparation

Meat (pork, beef, chicken und turkey) and sausage samples are grinded at room temperature with the Buchi Mixer B-400 before filling them to a Petri dish (see NIRFlex Solid package with Mixer and applications: N500-011 or N500-012).

The Ham application was done with a Fibre Optic Probe using whole pieces of Ham.

BÜCHI Labortechnik AG

CH - 9230 Flawil 1 T +41 71 394 63 63 F +41 71 394 65 65 buchi@buchi.com www.buchi.com

BÜCHI Labortechnik GmbH

D – 45127 Essen Freecall 0800 414 0 414 T +49 201 747 490 F +49 201 237 082 deutschland@buchi.com www.buchigmbh.de

BUCHI Sarl

F - 94656 Rungis Cedex T +33 1 56 70 62 50 F +33 1 46 86 00 31 france@buchi.com www.buchi.fr

BUCHI UK Ltd.

GB - Oldham OL9 9QL T +44 161 633 1000 F +44 161 633 1007 uk@buchi.com www.buchi.co.uk

BÜCHI Labortechnik GmbH

NL - 3342 GT Hendrik-Ido-Ambacht T +31 78 684 94 29 F +31 78 684 94 30 netherlands@buchi.com www.buchi.nl

BÜCHI Italia s.r.l.

I – 20090 Assago (MI) T +39 02 824 50 11 F +39 02 57 51 28 55 italia@buchi.com

BUCHI Hong Kong Ltd.

HK – Central T +852 2389 2772 F +852 2389 2774 china@buchi.com www.buchi.com.cn

BUCHI Shanghai

RC - 500052 Shanghai T +86 21 6280 3366 F +86 21 5230 8821 china@buchi.com www.buchi.com.cn

BUCHI (Thailand) Ltd.

T - Bangkok 10600 T +66 2 862 08 51 F +66 2 862 08 54 bacc@buchi.com www.buchi.com

Nihon BUCHI K.K.

J – Tokyo 110-0008 T +81 3 3821 4777 F +81 3 3821 4555 nihon@buchi.com www.nihon-buchi.jp

BUCHI India Private Ltd.

IND - Mumbai 400 055 T +91 22 667 18983 / 84 / 85 F +91 22 667 18986 india@buchi.com www.buchi.in

BUCHI Canada Ltd.

CDN-Thornhill, Ontario L4J6Z2 T +1 416 277 7407 F +1 905 764 5139 canada@buchi.com www.mybuchi.com

BUCHI Corporation

USA - New Castle, Delaware 19720 Toll Free: +1 877 692 8244 T +1 302 652 3000 F +1 302 652 8777 us-sales@buchi.com www.mybuchi.com

We are represented by more than 100 distribution partners worldwide. Find your local representative at

